


Instructions:

1. Connect the Cyclops to a computer using a mini-USB to USB cord.
2. Run the UM_Factory.exe file

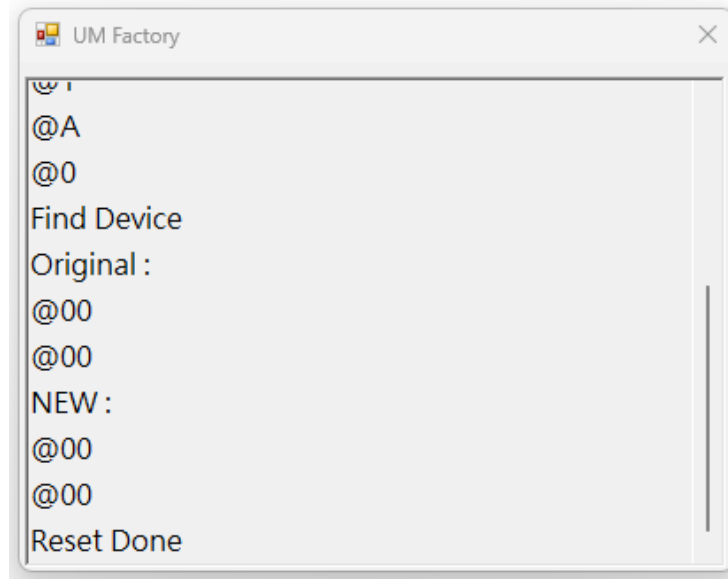

When you see **NEW:**, the process is complete.

@00
@00

After resetting the motor, please zoom in or out to achieve a field of view (FOV) of 8 mm and check the magnification displayed.

To confirm the correct magnification for the displayed FOV, you can calculate it manually using the following formula:

Monitor width (mm) / FOV (mm)

For example, with an FOV of 110 mm: 541 mm / 110 mm = 4.9x